

NEW SPACE FOR YOUR DATA-

Imaging Facility Disc Station

Imaging Facility Disc Station is a new space where Imaging Facility users can save and store their data from Spinning disc and Zeiss880 microscopes.

How to map the Disc Station from ZEISS880 microscope PC:

- Log in as LSM880\user to Windows (or as Asuch user)
- Select **Computer** from the Start menu
- Select **Map network drive**
- Name the drive as **I:**
- Select from history or type <\\ds-ueb-if.ueb.cas.cz/home>
- check checkbox → connect using different credentials
- Press **Finish** button
- log in as ASUCH user (your name and pswd)

The Imaging Facility Disc Station will is available as your new drive (I:) on microscope PC, you can save your data directly there.

- Before leaving disconnect the I: drive by right-mouse button on I: drive and select Disconnect

How to reach your data

- in ASUCH domain, map the Imaging Facility Disc Station on your PC analogically as above

From Windows computer out of ASUCH domain:

- type web address <http://ds-ueb-if.ueb.cas.cz>
- log in
- choose left top icon
- in File Station work with your files

From Mac:

- map disc using Finder, select on top **Go -> Connect to Server**. Fill in path *afp://ds-ueb-if.asuch.cas.cz/home* and log in as in ASUCH domain.

Assistance: Katka or Honza 1.435