

Zeiss Axio Imager + ApoTome2

HELP and training for new users:

Katka Malínská tel. 435 or 732 271 456

Start the system

- Turn on: **1) main switch**
2) fluorescence lamp HBO100 (if you will need it)
3) microscope (on the microscope body left)
4) PC as the last one
- User **CZC54304YV\zeiss** no password
- Start **Zen blue** and choose **ZEN Pro** mode

Acquisition

you can e.g. use saved configurations for individual fluorophores

Shift + F2 – live camera mode

F2 – capture image

Saving your data

I. Locally on D:\DATA USER\your name

II. or map imaging facility disc station:

<\\ds-ueb-if.ueb.cas.cz\home>

III. or map your P: drive disc station:

<\\ds.asuch.cas.cz\home>

(login as in asuch domain)

ApoTome structured illumination

- Insert ApoTome slider
- In Zen activate ApoTome function and choose number of grid positions, increase time exposition

Turn off the system

- Turn off:
- 1) PC**
 - 2) fluorescence lamp HBO100**
 - 3) microscope** (on the microscope body left)
 - 4) main switch**
 - 5) If you have used the ApoTome slider turn in out again**

Please let us know about any problems. Thank you.

Katka and Honza
tel: 435; 732271456